

# Microbiological and Particle Control of Compressed Gases

A review

FRANK PANOFEN, PhD DANIELE PANDOLFI



### Agenda

- Compressed Gases as an Essential Part of an EM Program
- Application overview
- Microbial Survival in Compressed Gases Under Fast Decompression to Normal Atmospheric Conditions
- Instrument Selection Criteria and ROI
- Conclusion



Without measurement there is no control

### Compressed Gases as an Essential Part of an EM Program





- Control of the environment in which pharmaceutical products are manufactured is a key element of Good Manufacturing Practices (GMP).
- Physical and microbial monitoring of manufacturing cleanrooms, Restricted-Access Barrier Systems (RABS), and isolators consists of clearly defined components.
- Absence of microbiological contamination is considered a critical quality attribute due to its potential to dramatically impact, directly or indirectly, the safety and/or the efficacy of the drug product.
- Compressed gas is considered a critical utility in the Pharma industry.
  - Direct product contact
  - Direct entry into clean environments



### Physical monitoring of manufacturing cleanrooms

### Components of an environmental monitoring program

Туре	Component	Characteristics	Purpose		
Physical	<ul> <li>Total particulate of room or enclosure air + Compressed Gases</li> </ul>	<ul> <li>Total inert and viable particles</li> <li>Active sampling</li> <li>Particle counter</li> <li>Decompression adaptor</li> </ul>	<ul> <li>Quality of environment</li> <li>Root-cause analysis for source of contamination</li> <li>Evaluation of compressed gas quality: Equal or better than air in the critical area</li> </ul>		
	<ul> <li>Differential pressures, positive pressure from a critical area to adjacent area</li> </ul>	<ul><li> 10 - 15 Pa difference</li><li> Sensor</li></ul>	<ul> <li>Prevents entry of contaminant particles from adjacent to critical clean area</li> </ul>		
	<ul><li>Air changes</li><li>Unidirectional air flow</li></ul>	• Sensor	Quality of environment		
	Temperature and     relative humidity	Thermometer,     hygrometer sensors	<ul> <li>Product quality and personnel comfort</li> </ul>		



### Microbial monitoring of manufacturing cleanrooms

### Components of an environmental monitoring program

Туре	Component	Characteristics	Purpose
Microbiological	<ul> <li>Viable counts of room or enclosure air &amp; compressed gases</li> </ul>	<ul> <li>Microbial counts by active (impactors and others) sampling</li> <li>Passive (settle plate) air sampling</li> <li>Sampling of gas by microbial gas sampling impactor with decompression adapter</li> </ul>	<ul> <li>Quality of environment</li> <li>Root-cause analysis for source of contamination</li> <li>Evaluation of compressed gas quality: Equal or better than air in the critical area</li> </ul>
inici o biological	<ul> <li>Viable counts on personnel</li> <li>Viable counts on surfaces</li> </ul>	<ul> <li>Effective personnel training and hygiene</li> <li>Microbial counts by contact plates and swabs</li> </ul>	<ul> <li>Contamination risk from operators</li> <li>Evaluation of aseptic techniques and practices</li> </ul>
		<ul><li>Microbial counts by</li><li>contact plates &amp; swabs</li></ul>	<ul> <li>Effectiveness of aseptic techniques and cleaning</li> <li>Disinfection of area and equipment</li> </ul>



Regulatory requirements

- There is a lack of specific guidelines/regulation for pharmaceutical applications.
- Key sources of information: FDA/GMP Guidelines
- Additional information may come from **ISO standard 8573**
- Key attributes for CG applications
  - Solid particles
  - Water
  - Oil
  - Microbiological burden





"A compressed gas should be of appropriate purity (e.g. from oil) and its microbiological and particle quality after filtration <u>should be equal or better than that of the air in the environment</u> <u>into which the gas is introduced</u>. Compressed gases such as air, nitrogen and carbon dioxide are often used in cleanrooms and are frequently employed in purging or overlaying."

(FDA aseptic filling guide)



- Part 1: Contaminants and purity classes
- Part 2: Test methods for aerosol oil content
- Part 3: Test methods for measurement of humidity
- Part 4: Test methods for solid particle content
- Part 5: Test methods for oil vapor and organic solvent content
- Part 6: Test methods for gaseous contaminant content
- Part 7: Test method for viable microbiological contaminant content
- Part 8: Test methods for solid particle content by mass concentration
- Part 9: Test methods for liquid water content





Solid particulate, water and oil concentration per ISO class										
		SOLID PARTICULATE			WATER		OIL			
ISO8573-1:2010 Class	Maximum number of particles per m3			Mass Concentration	Vapour Pressure Dewpoint	Liquid g/m³	Total Oil (aerosol liquid and vapour)			
	0.1 - 0.5 micron	0.5 - 1 micron	1 - 5 micron	mg/m <sup>3</sup>	°C		mg/m³			
0	As specified by the equipment user or supplier and more stringent than Class 1									
1	•NOTE: ISO 8573 classes are not identical to the 1 ISO 14644 classes.									
2										
3										
4										
5	<ul> <li>If you choose to use the 8573 concentration</li> </ul>									
6	limit table for your CG classification, make sure									
7	the appropriate particles sizes are considered,									
8										
9	respective to your ISO 14644 class limits.									
X	-	-	_				10			



Sampling point classification – Summary

- Typically, in an aseptic filling environment this table will not apply (equivalent to Class 0 in 8573-1).
- Manufacturers will need to define ther own specifications according EU/FDA GMP/ISO guidelines.
- In pharmaceutical production environments, the cleanroom user shall guarantee the compressed gas purity to meet or exceed the same cleanroom grade limits.
  - For example, a gas used in a Grade A/ISO 5 area should not exceed the Grade A/ISO 5 maximum permitted particle limit.



*ISO 8573-7 test method for viable microbiological contaminants* 

- ISO 8573 Part 7 provides a means of sampling, incubating and determining the number of microbiological particles.
  - Sampling method: Impaction (STA) together with the method given in ISO 8573-4
  - Isokinetic sampling of the air shall be carried out and reduced until it is within the range of the sampler as identified by the manufacturer.
  - Pressure reduction to atmospheric conditions and flow measurements shall be performed to establish compatibility with the manufacturer's recommendations or in accordance with ISO 8573-4.
  - Limits are not specified.


⇒Note: Investigations have shown that less than 50% of pharmaceutical plants WW use the impaction method (direct impaction or STA).



# Example #1: Microbial CG sampling setup

### MiniCapt Mobile CG Kit

- Special diffusion chamber mounted to sampler inlet
- Other end connected to the CG distribution lines
- Sampling at balanced flow
  - Compressed gas flow = sampling flow
- The impactor inlet pressure coincides with the pressure of the compressed gas kit diffusion chamber.
- Pressure regulator between the CG source & inlet
  - Safety
  - Regulate the flow rate of the gas into the diffusion chamber
- Solenoid valve connected downstream of the regulator allows synchronization of gas with the start of a sample.
- Available in 25, 50 or 100 LPM configurations
- Active monitoring of gas flow to detect changes

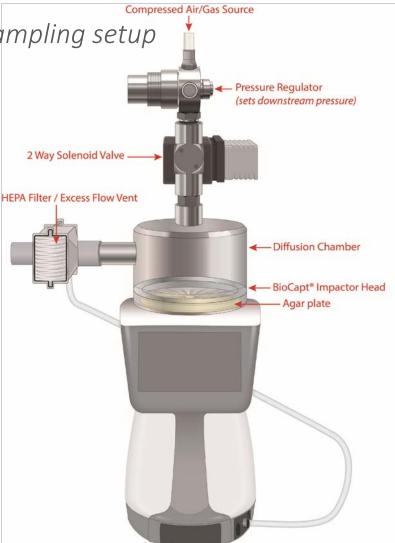



Example #2: Particle CG sampling setup

### Lasair III HPD – High Pressure Diffuser

- Provides a plug & play solution for particle contamination control in CG.
- No pressure adjustment required
- No additional calibration required
- The Critical Orifice principle guarantees a stable flow in accordance with the instrument flowrate requirement.
- The HEPA filter guarantees any contaminated exceeding gas pressure will not contaminate the cleanroom.



- Available in 28.3, 50 or 100 LPM configurations
- Accept gas pressure from 25 to 100 psi
- Designed for portable or fixed installation



### Frequency

- Initial qualification
  - 1 week daily
  - Monthly
- Routine monitoring
  - Quarterly
- Long term routine monitoring
  - Risk based approach
  - Annually to bi-annually

### Location

- Initial qualification
  - Every sample location
- Routine monitoring
  - Worst case locations based on initial qualification
- Long term routine monitoring
  - Risk based approach based on historical data



Without measurement there is no control

## Applications





Gas applications in aseptic environments used in pharmaceuticals & biotechnology

### **Example of Gas Applications in Pharmaceutical Industry**

#### Blanketing:

Maintain an inert atmosphere above a liquid or powdered product inside a container. Prevention of product degradation/corrosion, control of volatile emissions.

#### Aseptic Packaging:

Packaging of medical supplies

Maintain sterility and cleanliness of the product by purging oxygen from the packaging

#### **Pressure and Leak Testing:**

Inert gas to pressurize new, repaired, or modified tanks, pipelines, vessels, and process equipment in order to check their integrity and leak tightness.

#### **Pressure Transfer:**

Transfer liquid or powder products to and from containers without requiring pumps


#### Filtration, Separation and Purification

**Purging:** 

The use of inert gas to displace air, flammable vapors, and contaminants from containers



### Use of compressed gases in Pharmaceutical Industry % \*

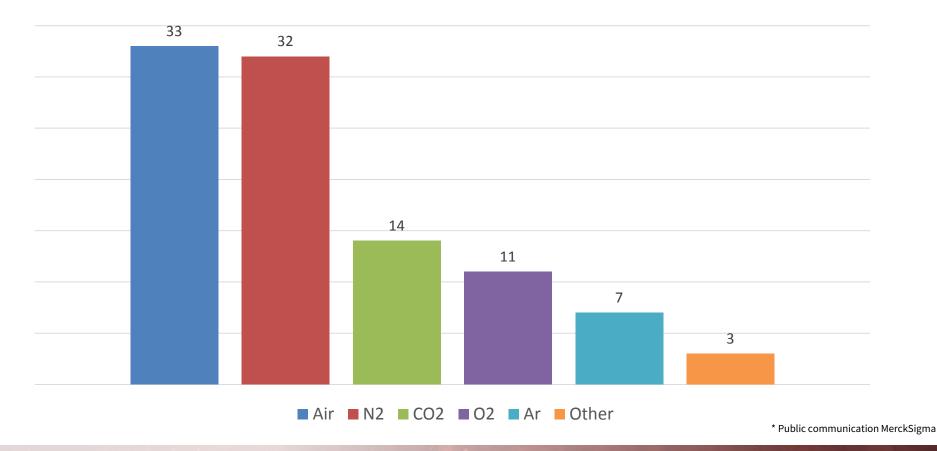

Confidential and proprietary

Without measurement there is no control



Without measurement there is no control

Microbial Survival in Compressed Gases Under Fast Decompression to Normal Atmospheric Conditions





Bacterial cells survive high pressures!

Bacterial cultures in food products require 2500- 3000 bar\* to inactivate<sup>3</sup> Lessened microbial growth & viability on metabolic processes in vivo are seen with pressures higher than 1000 bar\*<sup>5</sup>

Cells treated at pressures of <1000 bar\* showed **no significant loss** of viability<sup>3</sup>

Serratia and Carnobacterium spp Survive conditions similar to Mars (low O2, <0°C, 7 mbar)<sup>5</sup>

\*1 Bar = 0.1 Mpa = 14.5 PSI

- 3 FDA. (2014). Kinetics of Microbial Inactivation for Alternative Food Processing Technologies - High Pressure Processing (HPP). A report of the Institute of Food Technologists for the Food and Drug Administration of the U.S. Department of Health and Human Services. Updated December 2014.
- 5 Nicholson, W. (2013). Growth of Carnobacterium spp. from permafrost under low pressure, temperature, and anoxic atmosphere has implications for Earth microbes on Mars. pnas 110(2):666- 671.




### Results<sup>2</sup>

- 92% recovery using the MAS100 CG
- Bacterial cells maintained suitable growth, even up to 10 bar during the incubation.
- Compression cycle did not influence the viability of the tested microorganisms.
- ⇒Hypothesis from study: Sampling prior to decompression avoided potential microbial cell damage and lowered viability. This hypothesis is refuted by multiple studies.



*Microbial cells without gas vesicles easily survive fast decompression from 10 – 1 bar* 

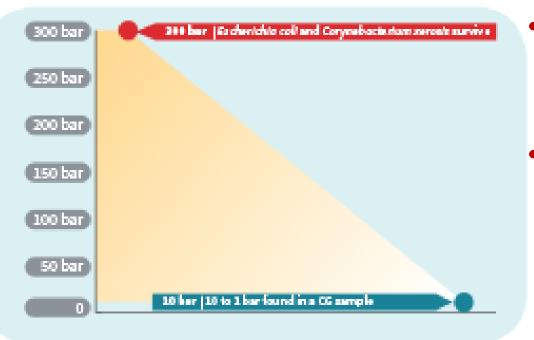


- Cells without gas vesicles were not harmed by decompression from up to 300 bar.<sup>4</sup>
- Escherichia coli and Corynebacterium xerosis survive rapid decompression from 300 bar (vs decompression from 10 to 1 bar found in a CG sampler]

4 Hemmingsen, B. B. & Hemmingsen, E. A. (1980). Rupture of the Cell Envelope by Induced Intracellular Gas Phase Expansion in Gas Vacuolate Bacteria. Journal of Bacteriology. 143 no. 2:841-846.



- Gram-negative, gas vacuolate bacteria (*M. aquaticus, P. pneumaticum, and M. glaucopis*) are known to be very susceptible to decompression.<sup>4</sup>
  - Populations of these cells were saturated with Ar, N<sub>2</sub> or He up to 100 bar. Gas phases of the vesicles remained intact.
  - Upon rapid decompression to atmospheric pressure, the vesicles expanded and ruptured the cells.
  - Viable counts indicated minimum pressures were between 25 50 bar.
  - Majority of the cell envelopes were ruptured at pressures between 50 - 100 bar

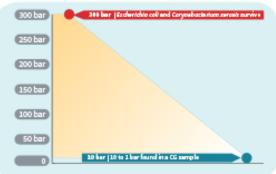


4 Hemmingsen, B. B. & Hemmingsen, E. A. (1980). Rupture of the Cell Envelope by Induced Intracellular Gas Phase Expansion in Gas Vacuolate Bacteria. Journal of Bacteriology. 143 no. 2:841-846.




Fast decompression - conclusion

• The analogy of bacterial cells and human divers is wrong!

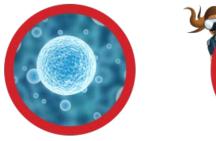



- Sensitive (vesicle containing) microorganisms are affected by rapid decompression.
- The viability is affected only when compression exceeds 25 bar to 1 bar.
- These conditions do not compare to realistic pharmaceutical settings.

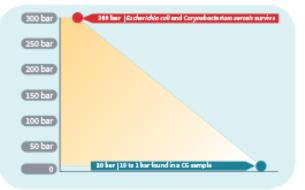



Without measurement there is no control

### Instrument Selection Criteria and ROI





MICROBIAL Instruments or adaptors from multiple vendors are available



MAS CG EX<sup>®</sup>, (MerckSigma/MBV)



RCS <sup>®</sup> CG Adaptor, (MerckSigma)



MiniCapt<sup>®</sup> Mobile, (PMS)



### Instruments or adaptors from multiple vendors are available



Lighthouse High Pressure Diffuser



PMS - High Pressure Diffuser III



MET ONE High Pressure Diffuser



Climet High Pressure Diffuser



Challenges of CG monitoring

- Compressed gases need to be tested directly on-line
  - Especially in critical areas
- Reality: Monitoring points are not always established close to the point of use
  - ⇒Size and weight of instrument mobile units
  - $\Rightarrow$ Ease of use
  - ⇒Software/data capabilities
  - ⇒Cleanroom protection: HEPA filtration



Low accessibility of sampling points



### Selection of Instruments - ROI

### Challenge

- Investment for rare measurement
- Full service program for a dedicated instrument (IQ/OQ + maintenance/calibration)

### Solution

- External service provider
- Multi-functional instrument
  - Only cost of accessory
  - Full service program
  - Routine monitoring
  - Flexibility




Without measurement there is no control

### Conclusions





### Conclusions

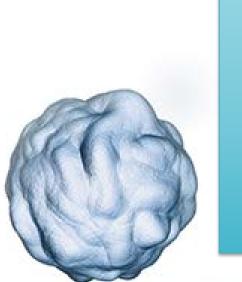

- Compressed gas monitoring is an essential part of an EM Program.
- Microorganisms survive hyperbaric or hydrostatic pressures from 1 to 10 bar.
- Decompression from 10 bar to normal atmospheric pressure does not harm microorganisms.
- Instrument selection and ROI calculations follow specific criteria of size/weight and flexibility of functionality.



# Thank you

# Frank Panofen fpanofen@pmeasuring.com

Daniele Pandolfi dpandolfi@pmeasuring.com

